

# LC/MS Application Note





# Determination of pharmaceuticals in water using online SPE coupled to liquid chromatography mass spectrometry

C. vom Eyser<sup>1</sup>, T. Teutenberg<sup>1</sup>, J. Tuerk<sup>1</sup>, Andreas Bruchmann<sup>2</sup>

<sup>1</sup> Institut für Energie- und Umwelttechnik e. V. IUTA (Institute of Energy and Environmental Technology) Bliersheimer Str. 58-60, 47229 Duisburg, Germany

<sup>2</sup> Axel Semrau GmbH & Co.KG, Stefansbecke 42, 45549 Sprockhoevel, Germany

Keywords: Pharmaceuticals, wastewater, online solid phase extraction, SPE, LC-MS

# **Overview Results**

- Monitoring of pharmaceuticals in surface and waste waters.
- conventional offline SPE LC-MS and online SPE LC-MS
- The challenging goal of 10 ng/L for monitoring was achieved by use of online SPE.
- Demonstration of a routine method for the online SPE LC-MS analysis of 24 pharmaceuticals in surface and wastewater.

#### Introduction

- The ubiquitous distribution of pharmaceuticals in different water bodies results in negative environmental effects. In Germany alone about 8'000 metric tons are drained into sewage systems every year. Roughly 90% of these compounds pass through treatment plants into different waterbodies.
- The low environmental concentrations require efficient enrichment and clean-up methods.
- Solid phase extraction (SPE) promises to achieve the monitoring target of 10 ng/L for each analyte.
- Automated enrichment and clean-up prior to liquid chromatography mass spectrometry (LC-MS) analysis using online SPE is demonstrated.

#### Methods



Figure 1: Instrument setup

## Offline SPE

Cartridge Phenomenex, Strata-X (polymeric reversed phase) Sample volume 100 mL - 1 L

Drying under nitrogen, Resolution in 1 mL water + 0.1% formic acid

#### Instrumentation online SPE

Sample prep	PAL RTC
Automated SPE	Axel Semrau CHRONECT µSPE Exchange Module
Cartridge type/format	HySphere C18 10 x 2 mm
LC	Agilent 1100
Column	Chromolith FastGradient Rp 18e 50-2 mm (Merck)
MS	Sciex API 3000 under Analyst 1.6.2
Control software	Chronos v. 4.4

#### Results

- C18 material was the best compromise between enrichment capability and peak shape for the analysis of the pharmaceuticals analyzed here.
- Enrichment factors of up to 200 (equal to 10 mL injection) resulted in good recovery rates.
- The recovery for conventional offline SPE and online SPE was comparable. The recovery of most analytes is between 80 and 120% (fig. 2)



Fig.2: Relative recovery and reproducibility for the 24 compounds analyzed.

- The average reproducibility for online SPE (5.1%) is better than for offline SPE (27%, see error bars in fig. 2)
- The analysis time/sample was minimized by efficient overlapping of tasks with an intelligent software algorithm (Chronos)



Fig.3: Comparison of time consumption for processing of 3 samples using offline SPE, online SPE and staggered online SPE.

• Applying online SPE the limit of detection (LOD) improved for example from 270 ng/mL to 3 ng/mL for metoprolol (fig. 4)



(Data from Sciex API 3000.)

Fig.4: A: Extracted ion chromatogram of a 10 mL online SPE sample of metoprolol (0.025 ng/mL) B: Extracted ion chromatogram of a 50 μL direct injection sample of metoprolol (5 ng/mL)

- The limit of detection is improved by a factor of 66 on average by enrichment via online SPE.
- For most pharmaceuticals the LOD is below the monitoring target of 10 ng/L (fig. 4) even though an older MS system was used.



Fig.5: Comparison of the limits of detection (LOD) for the compounds analyzed between online SPE (green) and direct injection (blue). For 22 of the 24 compounds the desired LOD < 10 ng/L was achieved with online SPE.

## Conclusions

- Online SPE is a suitable routine tool for the automated enrichment of pharmaceuticals from wastewater.
- Recovery and reproducibility of online SPE are comparable to or better than conventional offline SPE.
- The monitoring target of 10 ng/L is achieved for 22 of 24 pharmaceuticals, even with an older MS system.
- Chronos as control software allows to overlap SPE process and MS-analysis, resulting in a runtime of only 162 min for 10 samples.
- The automated process improves process safety and minimizes labor cost.

#### Legal Statements

CTC Analytics AG reserves the right to make improvements and/or changes to the product(s) described in this document at any time without prior notice.

CTC Analytics AG makes no warranty of any kind pertaining to this product, including but not limited to implied warranties of merchantability and suitability for a particular purpose.

Under no circumstances shall CTC Analytics AG be held liable for any coincidental damage or damages arising as a consequence of or from the use of this document.

© 2016 CTC Analytics AG. All rights reserved. Neither this publication nor any part hereof may be copied, photocopied, reproduced, translated, distributed or reduced to electronic medium or machine readable form without the prior written permission from CTC Analytics AG, except as permitted under copyright laws.

CTC Analytics AG acknowledges all trade names and trademarks used as the property of their respective owners.

PAL is a registered trademark of CTC Analytics AG | Switzerland 6 | IngeniousNews 06/2016

#### Imprint

Date of print: 12.2016

CTC Analytics AG Industriestrasse 20 CH-4222 Zwingen Switzerland T +41 61 765 81 00 Contact: info@ctc.ch